


### For PMA.core :

Host:  Username:  Password:  Folder:  

status: Trying to connect...  
status: Resolving address of   
status: Connecting to   
status: PMA.core v2.0.0.1686  
status: Connection established...  
status: Loading folder on PMA.core...  
status: Folder successfully loaded...

- Host : PMA.core instance to connect to, it should be in this form “http(s)://xxxxx/pma.core.2/”
- Username & Password : connection credentials
- Folder : Remote directory on server to navigate to if connection is successful

### For My Pathomation :

Email:  Password:  Folder:  

status: Trying to connect to the cloud...  
status: Resolving address of   
status: Connecting to   
status: PMA.core v2.0.0.1696  
status: Connection established...  
status: Loading user's folder on the cloud...  
status: Folder successfully loaded...

- Email : The email address used to subscribe to Pathomation cloud.
- Password : Password received on inbox once the subscription & payment are validated.
- Folder : disabled text field, end user can only connect to its own subscription folder.

From:  
<https://docs.pathomation.com/PMA.transfer/> - **PMA.transfer**

Permanent link:  
[https://docs.pathomation.com/PMA.transfer/doku.php?id=connection\\_bar\\_logging\\_text\\_area&rev=1589528638](https://docs.pathomation.com/PMA.transfer/doku.php?id=connection_bar_logging_text_area&rev=1589528638)

Last update: **2020/05/16 05:14**

