


	Slidename	Slidesize	FileType	Last Modified
↑	..			
📁	brightfield		File folder	
📁	testtempwim2		File folder	
📁	WG26		File folder	
🖼️	CMU-1.mrxs	565,102,675	MRXS File	Sun Dec 15 11:46:31 GMT+01:00 2019
🖼️	CMU-1.svs	177,552,579	SVS File	Sun Dec 15 11:47:06 GMT+01:00 2019

Initiated via mouse right click within the table view borders, it's composed of three menu items : "Rename folder", "Delete folder", "Download folder".

Rename folder :

Click on this option pops up the following window :

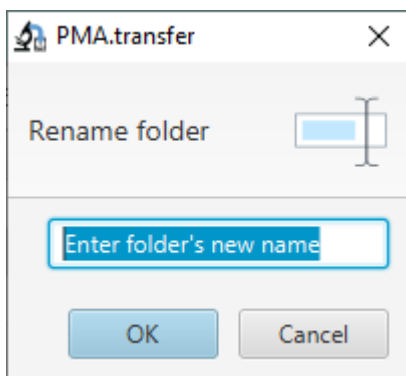

It's only permitted to rename empty folders and the new name shouldn't be already used on a neighboring folder.

Delete folder :

It's only permitted to rename an empty folder. If it's the case, a confirmation window asks for delete confirmation, once confirmed the folder is removed completely from the server (warning : this action cannot be undone).

Download folder :

Has been discussed in details earlier on this tutorial.

From: <https://docs.pathomation.com/PMA.transfer/> - PMA.transfer

Permanent link: [https://docs.pathomation.com/PMA.transfer/doku.php?id=context\\_menu\\_for\\_folders\\_remote&rev=1589592873](https://docs.pathomation.com/PMA.transfer/doku.php?id=context_menu_for_folders_remote&rev=1589592873)

Last update: 2020/05/16 05:14

