

There are two options to launch a download :

- Select a slide on local site (Table view to the left) then click on menu “Transfer” on the toolbar then select menu item “Download slide”
- Mouse right click inside the left table view borders, then select “Download slide to local folder” context menu item.

Download of a folder (recursively) :

The download of a folder can be performed via a click on context menu item “download folder to local site” . Once successfully initiated, The folder and its sub-directories are sub-consequently created on local side then the download of contained slides is operated to corresponding destinations.

From:

<https://docs.pathomation.com/PMA.transfer/> - **PMA.transfer**

Permanent link:

[https://docs.pathomation.com/PMA.transfer/doku.php?id=initiating\\_a\\_download&rev=1589595289](https://docs.pathomation.com/PMA.transfer/doku.php?id=initiating_a_download&rev=1589595289)

Last update: **2020/05/16 05:14**

