


# Menu bar

The menu bar is composed of five menu items : “File”, “Edit”, “View”, “Transfer”, “Help”




## File

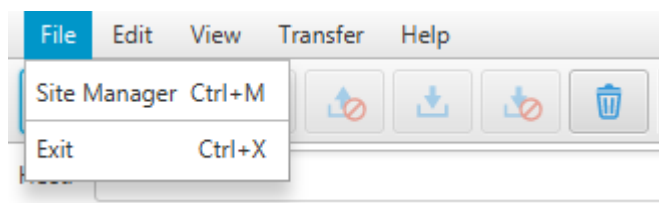


- Click on “Site Manager” launches Site manager window
- Click on “Exit”, shuts down PMA.Transfer

## Edit

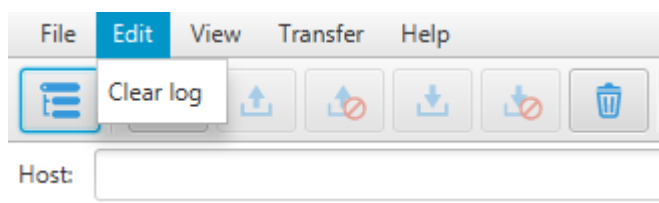


- Click on “Clear log”, clears logging messages in the text area below connection text fields

## View

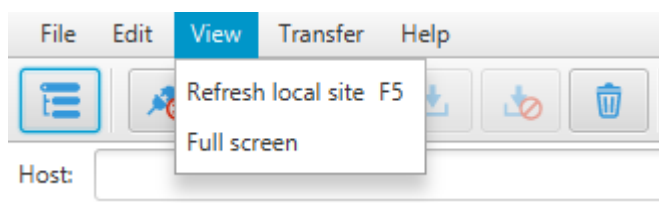


- Click on “Refresh local site”, refreshes the table view to the left related to content on end user’s local computer.
- Click on “Full screen” switches to full screen display.

## Transfer

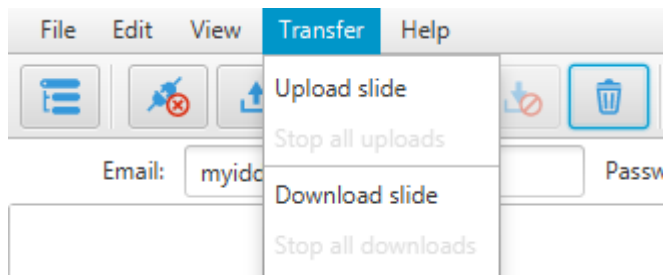

- Click on “upload slide” : Enabled when a slide on local HDD is selected, it initiates an upload to the remote site PMA.transfer is actually connected to (a couple of checks are operated before an upload can be started : duplicate slide name, slide already being uploaded, no connection...)
- Click on “Stop all uploads” : Enabled when there are active uploads running, it aborts all current running uploads.
- Click on “Download slide” : Enabled when a slide on remote site is selected, it initiates a download to the selected destination on local site (a couple of checks are operated before a download can be started : creation of destination folders, duplicate slide name, slide already being downloaded, no connection...)
- Click on “Stop all downloads” : Enabled when there are active downloads running, it aborts all current running downloads.


## Help

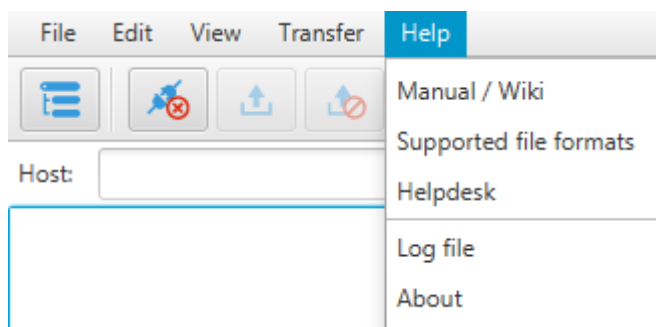

- Click on “Manual” downloads latest version of PMA.transfer manual.
- Click on “Supported file formats” loads a table view of the supported file formats by PMA.transfer.

Supported file formats					
vendor	extension	compression	brightfield	fluoro	zstack
TIFF	.tif, .tiff	JPEG, JPEG2000, LZW, Deflate, Raw, RLE	✓		
JPEG	.jpeg, .jpg	JPEG	✓		
JPEG 2000	.jp2	JPEG2000	✓		
PNG	.png	PNG	✓		
Olympus VSI	.vsi	lossless JPEG, JPEG, Raw	✓	✓	
Olympus WebView	.xml	JPEG	✓		
Ventana / Roche BIF	.bif	JPEG	✓		✓
Hamamatsu VMS	.vms	JPEG	✓		
Hamamatsu NDPI	.ndpi	JPEG	✓	✓	✓
Hamamatsu NDPIS (Image Set)	.ndpis	JPEG	✓	✓	✓
Hamamatsu DICOM	.dcm	JPEG, JPEG2000	✓		
Huron Technologies	.tif	JPEG	✓		
3DHistech MRXS	.mrxs	JPEG, PNG, BMP	✓	✓	✓
Aperio / Leica SVS	.svs	JPEG, JPEG2000	✓	✓	✓
Aperio / Leica AFI (Image Set)	.afi	JPEG, JPEG2000	✓	✓	✓
Aperio / Leica CWS	.cws	JPEG	✓		

- Click on “Helpdesk” redirects to a web interface to create a support ticket.


Guest User | [Sign In](#)

[Support Center Home](#)
[Open a New Ticket](#)
[Check Ticket Status](#)

## Welcome to the Support Center

In order to streamline support requests and better serve you, we utilize a support ticket system. Every support request is assigned a unique ticket number which you can use to track the progress and responses online. For your reference we provide complete archives and history of all your support requests. A valid email address is required to submit a ticket.

[Open a New Ticket](#)
[Check Ticket Status](#)

- Click on “Log file” opens up the current log file in notepad in addition to log files's location
- Click on “About” pops up an interface regarding general information about PMA.transfer



From:

<https://docs.pathomation.com/pma.transfer/> - **PMA.transfer**

Permanent link:

[https://docs.pathomation.com/pma.transfer/doku.php?id=menu\\_bar](https://docs.pathomation.com/pma.transfer/doku.php?id=menu_bar)

Last update: **2021/06/18 17:09**

