



- Click on “upload slide” : Enabled when a slide on local HDD is selected, it initiates an upload to the remote site PMA.transfer is actually connected to (a couple of checks are operated before an upload can be started : duplicate slide name, slide already being uploaded, no connection...)
- Click on “Stop all uploads” : Enabled when there are active uploads running, it aborts all current running uploads.
- Click on “Download slide” : Enabled when a slide on remote site is selected, it initiates a download to the selected destination on local site (a couple of checks are operated before a download can be started : creation of destination folders, duplicate slide name, slide already being downloaded, no connection...)
- Click on “Stop all downloads” : Enabled when there are active downloads running, it aborts all current running downloads.

From:
<https://docs.pathomation.com/PMA.transfer/> - **PMA.transfer**

Permanent link:
<https://docs.pathomation.com/PMA.transfer/doku.php?id=transfer&rev=1589528002>

Last update: **2020/05/16 05:14**

