

# System Logs

In the System Logs page of the System section of the administrators toolbar, you can see a list of errors that have been triggered.

Here you can see:

- 1. Host name
- 2. Error code
- 3. Type
- 4. Error description
- 5. User that triggered the error
- 6. Date of error
- 7. Time of error

The number of logs per page can be adjusted.

Error Log for pma.control.2 on

[RSS FEED](#) | [RSS DIRECT](#) | [DOWNLOAD LOG](#) | [HELP](#) | [ABOUT](#)

Errors 1 to 15 of total 969 (page 1 of 65). Start with 10, 15, 20, 25, 30, 50 or 100 errors per page.

Host	Code	Type	Error	User	Date	Time
	404	Http	Exception of type 'System.Web.HttpException' was thrown. <a href="#">Details...</a>	pma_admin	1/3/2024	9:54 AM
	404	Http	Exception of type 'System.Web.HttpException' was thrown. <a href="#">Details...</a>	User	12/22/2023	4:50 PM
	0	Web	The remote server returned an error: (500) Internal Server Error. <a href="#">Details...</a>		12/14/2023	12:47 PM
	404	Http	Exception of type 'System.Web.HttpException' was thrown. <a href="#">Details...</a>	b_user1	10/31/2023	2:34 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:55 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:55 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:55 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:55 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:55 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:55 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:54 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:52 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:52 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:34 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:31 PM
	404	Http	The controller for path  was not found or does not implement IController. <a href="#">Details...</a>	pma_admin	10/23/2023	1:30 PM

[Next errors](#)

Powered by [ELMAH](#), version 1.2.14705.955. Copyright (c) 2004, Atif Aziz. All rights reserved. Licensed under [Apache License, Version 2.0](#). Server date is Wednesday, 03 January 2024. Server time is 15:00:20. All dates and times displayed are in the W. Europe Standard Time zone. This log is provided by the XML File-Based Error Log.

From:  
<https://docs.pathomation.com/pathotrainer/> - **Pathotrainer**

Permanent link:  
[https://docs.pathomation.com/pathotrainer/doku.php?id=system\\_logs](https://docs.pathomation.com/pathotrainer/doku.php?id=system_logs)

Last update: **2024/01/03 17:21**