


Annotation Import and Export and 3rd Party Annotations

PIMS LS allows you to import and export annotations in a variety of formats, including Well-known text (the format used by PMA.core internally), Visiopharm mld files, Indica Labs HALO annotations files, Definiens/Aperio xml files, and 'plain' csv files.

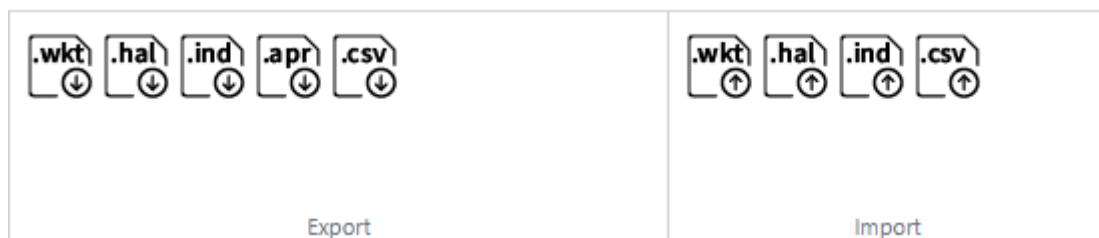


Export

You can export slide annotations in different formats including Well-known text (the format used by PMA.core internally), Visiopharm mld files, Indica Labs HALO annotations files, Definiens/Aperio xml files, and 'plain' csv files.

Import

User can also import/upload annotations through PIMS LS. Once you click on the import button for the format you want to import, a menu will appear prompting you to select the file that contains annotation data.

Once a valid file is uploaded annotations are added to the slide.

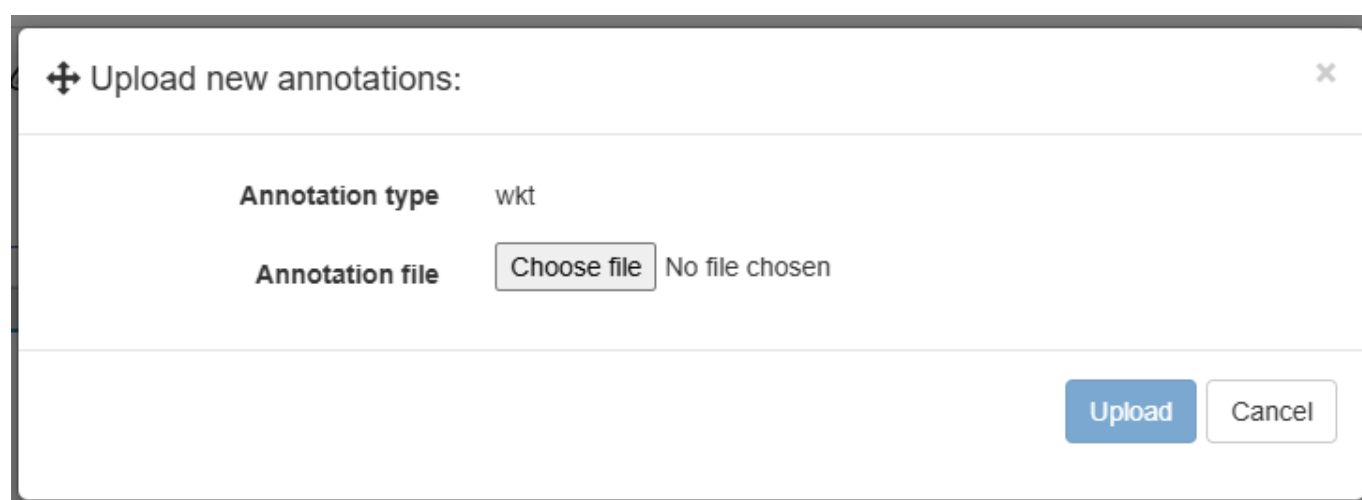

Last
update: 2024/11/29 11:38 annotations_import_export https://docs.pathomation.com/pimsls/3.0.0/doku.php?id=annotations_import_export&rev=1732869525

From:
<https://docs.pathomation.com/pimsls/3.0.0/> - **PIMS LS 3.0.0**

Permanent link:
https://docs.pathomation.com/pimsls/3.0.0/doku.php?id=annotations_import_export&rev=1732869525

Last update: **2024/11/29 11:38**

