


Version retrieval (automation)

For monitoring purposes, it is useful to determine the PMA.studio version used through a back-end API call. All of [Pathomation's platform](#) components have this option, actually.

In order to determine the installed version of PMA.studio, you can issue the call:

<https://server/pma.studio/api/json/getversioninfo>

The result is a JSON string formatted as A.B.C.D, with A, B, C being numeric. D can be either numeric or alphanumeric.

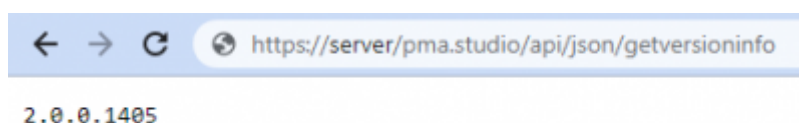

Reporting this version number to Pathomation when [asking for support through our helpdesk](#) is always helpful, too.

From:

<https://docs.pathomation.com/pimsls/3.0.0/> - **PIMS LS 3.0.0**

Permanent link:

https://docs.pathomation.com/pimsls/3.0.0/doku.php?id=auto_version

Last update: **2022/01/25 14:05**

