


# Home: Folder Tools Group


The Folder Tools group in the Home tab contains controls for searching & navigating folders

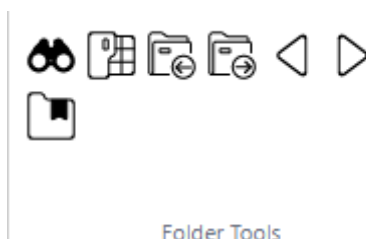

The icons, from left to right

## Top Row

1. - Search icon - [read more about searching here](#)
2. - Open current folder in grid - opens the menu below:

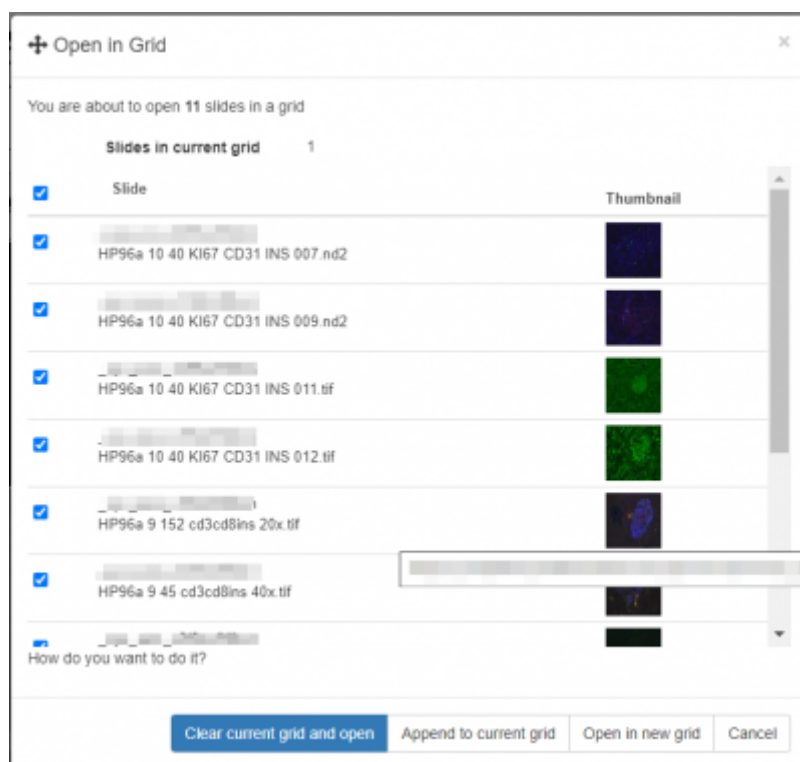

Here you can choose images from the currently highlighted folder in the slide browser to open into your current grid, or make a new grid.

3. - Previous folder - selects and opens the previous folder
4. - Next folder - selects and opens the next folder
5. - Previous file - opens previous image in folder
6. - Next file - opens next image in folder

From:

<https://docs.pathomation.com/pimsls/3.0.1/> - **PIMS LS 3.0.1**

Permanent link:

[https://docs.pathomation.com/pimsls/3.0.1/doku.php?id=folder\\_tools&rev=1731586987](https://docs.pathomation.com/pimsls/3.0.1/doku.php?id=folder_tools&rev=1731586987)

Last update: **2024/11/14 15:23**

