


Home: Grid Tool

The Grid tool contains icons for synchronising different viewports in the grid, clearing the grid, and opening new grid tabs.

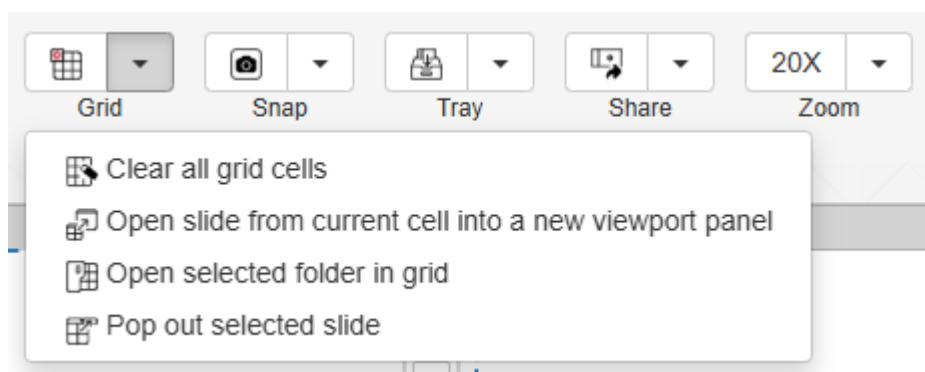

- Clear all grid cells - clears all images from current grid
- Opens current folder in grid - opens the menu below
- Opens slide from current cell into a new viewport panel
- Pops out selected slide in a new window.
- Clear current cell - clears image from currently selected grid cell

From:

<https://docs.pathomation.com/pimsls/3.0.1/> - **PIMS LS 3.0.1**

Permanent link:

<https://docs.pathomation.com/pimsls/3.0.1/doku.php?id=grid&rev=1746901175>

Last update: **2025/05/10 21:19**

