

## Available ribbon commands

See also the [tab inventory page](#).

Ribbon section	Command name	Parameters
File menu	custom-panel-algo-ichi	props=
	custom-panel-algo-ni	props=
	custom-panel-ai-prov-mindpeak	props=
	custom-panel-ai-prov-spintellx	props=
	file	
	about-us	
	show-eula	
	manual	
	sign-out	
Ribbon section	Command name	Parameters
Home tab	simple-layout	
	detailed-layout	
	detailed-layout	
	viewport-toggle-overview	
	viewport-toggle-slide-path	
	viewport-toggle-barcode	
	viewport-toggle-rotation-flip	
	add-tab	
	reset-layout	
	viewport-toggle-zoom	
	viewport-toggle-scalebar	
	viewport-toggle-label	
	panel-toggle-color	
	panel-toggle-search	
	open-folder-in-grid	
	prev-directory	
	next-directory	
	prev-image	
	next-image	
	add-folder-bookmark	
	reset-grid	
	clear-cell	
	open-in-tab	
	popout-info	
	snapshot	
	quick-snapshot	
	toggle-toolbar	
	increase-zoom	
	decrease-zoom	
	custom-zoom-1x	pma-magnification=
custom-zoom-2x	pma-magnification=	

<b>Ribbon section</b>	<b>Command name</b>	<b>Parameters</b>
	<a href="#">custom-zoom-5x</a>	pma-magnification=
	<a href="#">custom-zoom-10x</a>	pma-magnification=
	<a href="#">custom-zoom-20x</a>	pma-magnification=
	<a href="#">custom-zoom-40x</a>	pma-magnification=
	<a href="#">panel-toggle-color</a>	
	<a href="#">send-slide-to-tray</a>	
	<a href="#">send-to-tray</a>	
	<a href="#">send-dir-to-tray</a>	
	<a href="#">panel-toggle-tray</a>	
	<a href="#">share-slide-tray</a>	
	<a href="#">quick-share-folder</a>	
	<a href="#">quick-share-slide</a>	
	<a href="#">quick-share-slide-roi</a>	
	<a href="#">quick-share-grid</a>	
	<a href="#">quick-share-grid-rois</a>	
	<a href="#">quick-share-thumbnail</a>	
	<a href="#">share-slide</a>	
<b>Ribbon section</b>	<b>Command name</b>	<b>Parameters</b>
<a href="#">Grid tab</a>	<a href="#">toggle-1x1-grid</a>	
	<a href="#">toggle-2x2-grid</a>	
	<a href="#">toggle-3x3-grid</a>	
	<a href="#">sync-views-toggle</a>	
	<a href="#">reset-grid</a>	
	<a href="#">clear-cell</a>	
	<a href="#">open-in-tab</a>	
	<a href="#">viewport-add-row</a>	
	<a href="#">viewport-remove-row</a>	
	<a href="#">viewport-remove-row</a>	
	<a href="#">viewport-remove-column</a>	
	<a href="#">flip-rotate-menu</a>	
	<a href="#">reset-flip</a>	
	<a href="#">flip-horizontally</a>	
	<a href="#">flip-vertically</a>	
	<a href="#">reset-rotation</a>	
	<a href="#">rotate-clockwise-90</a>	
	<a href="#">rotate-counter-clockwise-90</a>	
	<a href="#">rotate-180</a>	
	<a href="#">zoom-in</a>	
	<a href="#">panel-toggle-annotations</a>	
	<a href="#">save-annotations</a>	
	<a href="#">toggle-annotations</a>	
	<a href="#">toggle-annotation-text</a>	
	<a href="#">toggle-fingerprint</a>	
<b>Ribbon section</b>	<b>Command name</b>	<b>Parameters</b>
<a href="#">Annotations tab</a>	<a href="#">Cursor</a>	

<b>Ribbon section</b>	<b>Command name</b>	<b>Parameters</b>
	delete-annotation	
	delete-all-annotations	
	edit-annotation-text	
	merge-selected-annotations	
	union-selected-annotations	
	modify-selected-annotation	
	Point	
	Circle	
	Line	
	Arrow	
	Ellipse	
	Rectangle	
	Landmark1	
	MultiPoint	
	Polygon	
	ClosedFreehand	
	CompoundFreehand	
	Freehand	
	Wand	
	Brush	
	export-to-wkt	
	export-to-mld	
	export-to-annotations	
	export-to-aperio	
	export-to-csv	
	import-from-wkt	
	import-from-mld	
	import-from-halo	
	import-from-csv	
<b>Ribbon section</b>	<b>Command name</b>	<b>Parameters</b>
Pre-set tab	preset-Circle0237mm2	pma-shape, pma-size-a, pma-size-b
	preset-Circle1mm	pma-shape, pma-size-a, pma-size-b
	preset-Circle2mm	pma-shape, pma-size-a, pma-size-b
	preset-Circle1cm	pma-shape, pma-size-a, pma-size-b
	preset-Square5mm	pma-shape, pma-size-a, pma-size-b
	preset-Square1cm	pma-shape, pma-size-a, pma-size-b
	preset-Tumor	pma-classification, pma-shape, pma-color
	preset-Macro	pma-classification, pma-shape, pma-color
<b>Ribbon section</b>	<b>Command name</b>	<b>Parameters</b>
Conferencing tab	create-conference	
	join-conference	
	start-conference	
	pause-conference	
	stop-conference	
	toggle-everyone-control-conference	

Ribbon section	Command name	Parameters
	<a href="#">broadcast-conference</a>	
	<a href="#">panel-toggle-conference</a>	
	<a href="#">panel-toggle-forms</a>	
	<a href="#">panel-toggle-links</a>	
	<a href="#">panel-toggle-navigation</a>	
	<a href="#">panel-toggle-bookmarks</a>	

From:

<https://docs.pathomation.com/pimsls/3.0.1/> - **PIMS LS 3.0.1**

Permanent link:

[https://docs.pathomation.com/pimsls/3.0.1/doku.php?id=ribbon\\_commands&rev=1680715695](https://docs.pathomation.com/pimsls/3.0.1/doku.php?id=ribbon_commands&rev=1680715695)

Last update: **2023/04/05 20:28**

