


# Ribbon layout (administration)

This admin page allows you to setup default Ribbon layout that consists of different tabs and controls. Each tab consists of different controls associated to it and you can configure appearance of tabs and controls using this page.

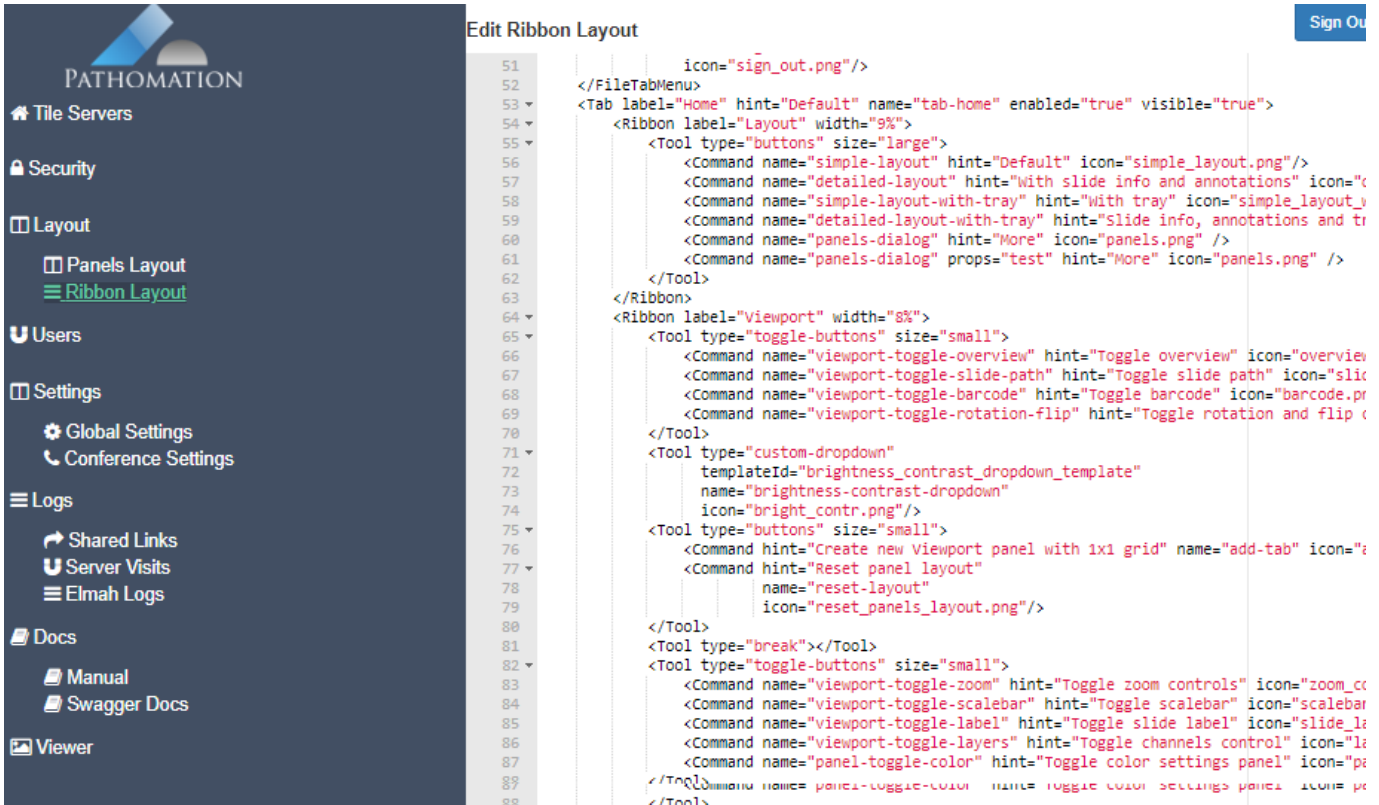

The screenshot shows the 'Edit Ribbon Layout' page in the Pathomation administration interface. The left sidebar contains a navigation menu with categories like Tile Servers, Security, Layout, Users, Settings, Logs, Shared Links, Server Visits, Elmah Logs, Docs, and Viewer. The 'Layout' section is expanded, showing 'Panels Layout' and 'Ribbon Layout'. The main content area displays the 'Edit Ribbon Layout' page with a code editor showing XML configuration for the ribbon layout. The code defines a 'Home' tab with various tools and commands, including 'viewport-toggle-rotation-flip', 'viewport-toggle-zoom', and 'panel-toggle-color'.

An [exhaustive list of all available commands](#) is available elsewhere on this site.

Below are some of the tabs and corresponding controls used in PMA.Studio.

**Home tab** has 8 sections:

1. View
2. Grid
3. Snap
4. Magnify
5. Directory
6. Slide
7. Share
8. Search



**Grid tab** supplies instruments to manage grid and it has 6 sections:

1. ItemSize
2. Synchronize
3. Rows/Columns
4. Slide

- 5. Snapshot
- 6. Share



**Annotations tab (available in PMA.Studio)** supplies instruments for working with annotations and has 4 sections:

- 1. Annotations
- 2. Shapes
- 3. Export
- 4. Import




**Pre-set Annotations tab (available in PMA.Studio)** consists of 3 sections:

- 1. Annotations
- 2. Stamps
- 3. Protocols




**Conferencing tab** has 2 sections:

- 1. Conferencing
- 2. Server

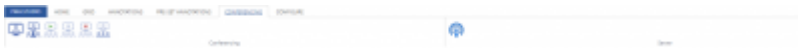

**Configure tab** has 4 sections:

- 1. Server
- 2. Panels
- 3. Pre-set layouts
- 4. Viewport

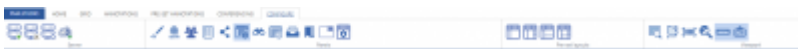

From:  
<https://docs.pathomation.com/pimsls/3.1.0/> - **PIMS LS 3.1.0**

Permanent link:  
[https://docs.pathomation.com/pimsls/3.1.0/doku.php?id=admin\\_ribbon&rev=1674470758](https://docs.pathomation.com/pimsls/3.1.0/doku.php?id=admin_ribbon&rev=1674470758)

Last update: **2023/01/23 13:45**

