


# Other Tools

Home tab also consists of independent icons like, Synx, grid overlays, toggle colors or search panel tools.

## Synchronize Tool

The Synchronize tool contains tools for turning on synchronized viewing and some grid management tools.

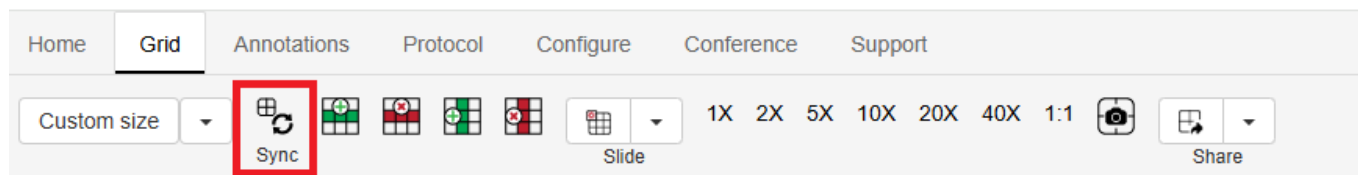



## Synchronized viewing

Turning on the Sync icon when you have 2 or more cells open in your grid will trigger synchronized viewing, you can see a quick video of this below:

## Grid Overlay Tools

Grid overlay tool allows user to add overlay to the opened slide. It can be of different sizes and colors. There are icons for both adding or removing overlay.

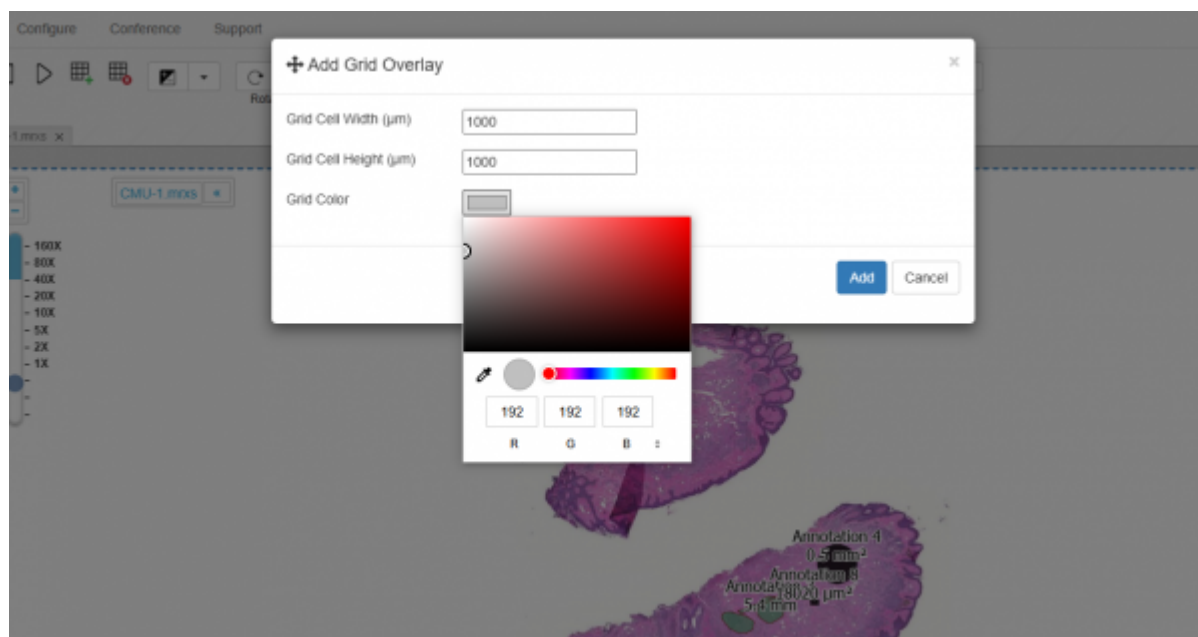

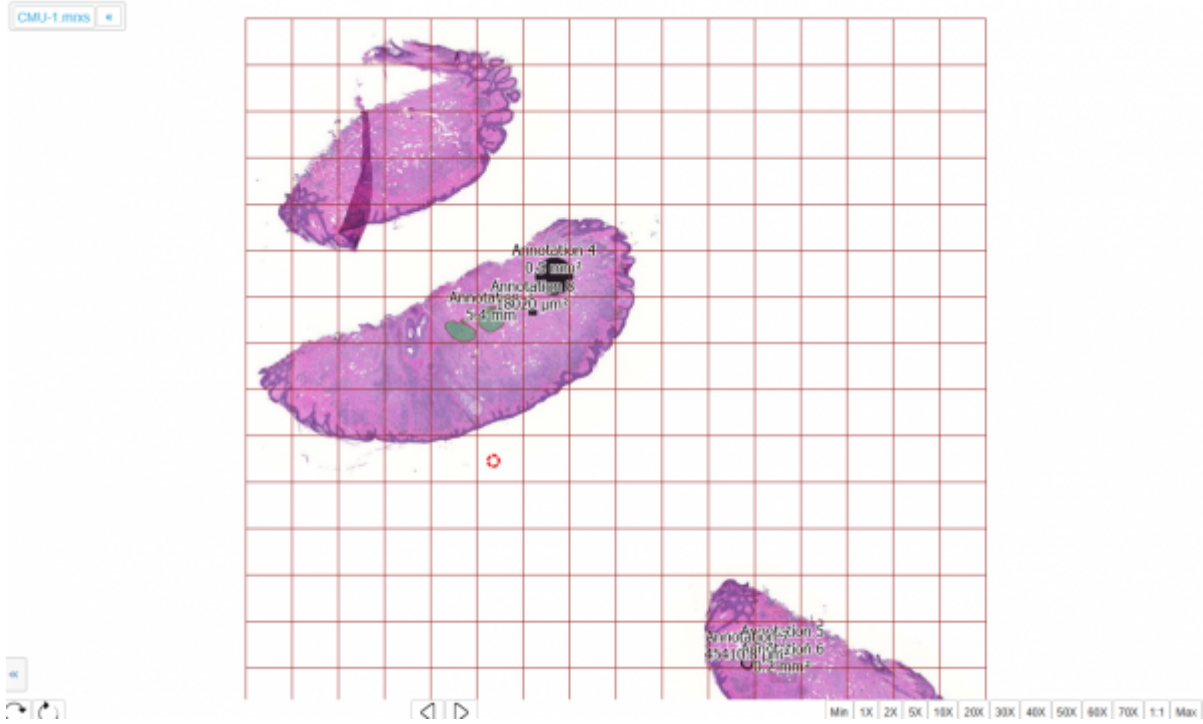

### Toggle panels

- Toggle color setting panel - [for more details](#)
- Search panel - [for more details](#)

From:  
<https://docs.pathomation.com/pimsls/3.1.0/> - **PIMS LS 3.1.0**

Permanent link:  
[https://docs.pathomation.com/pimsls/3.1.0/doku.php?id=other\\_icons&rev=1747061759](https://docs.pathomation.com/pimsls/3.1.0/doku.php?id=other_icons&rev=1747061759)

Last update: **2025/05/12 17:55**

