

The Forms Panel

The Forms panel is the place to input and review additional data about your slides

You can view a quick video guide on how to use the forms panel below:

[pimsls_formshowto.mp4](#)

Forms comprise of question/fields that you define.

Forms are completely customizable with many different datafields including:

- Numerical
- Freetext
- Checkbox
- Multiple choice
- Dropdown menu
- Percentage
- Date

Forms can be assigned at both the slide level, and at the folder / directory level so that you can have case-level and project-level data.

Please note forms need to be defined / created in PMA.core. Please ask your administrator if you don't have the form(s) you need.

From:

<https://docs.pathomation.com/pimsls/3.1.1/> - **PIMS LS 3.1.1**

Permanent link:

https://docs.pathomation.com/pimsls/3.1.1/doku.php?id=panel_forms

Last update: **2025/01/10 12:10**

