


Instances

This page shows a list of configured instances of PMA.core. An instance of PMA.core is an installation of PMA.core that share the same database as part of a group of Instances. This section is particular relevant in the context of [geo-replication scenarios](#).

Usually each instance is a replica of PMA.core at a different server and location. You can set an instance as the current running in this page by clicking on the *Edit → Set current* button

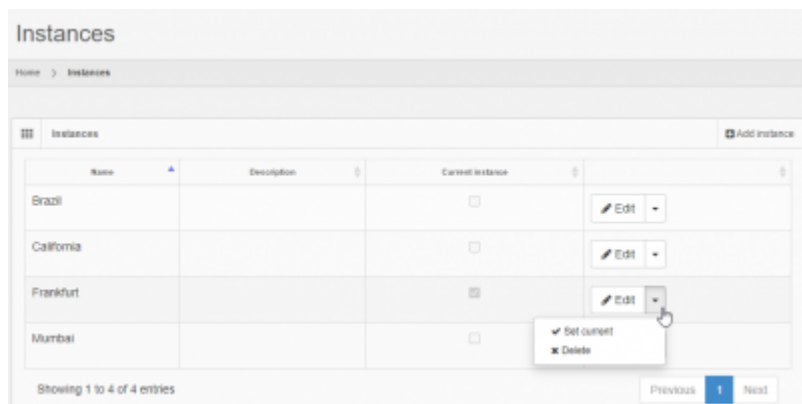


Name	Description	Current instance	Edit
Brazil		<input type="checkbox"/>	Edit
California		<input type="checkbox"/>	Edit
Frankfurt		<input checked="" type="checkbox"/>	Edit
Mumbai		<input type="checkbox"/>	Edit

Showing 1 to 4 of 4 entries

Previous 1 Next

Adding a new instance

Clicking on the *Add new instance* button will take you to the page to add a new Instance of PMA.core. The only needed information about an instance is a unique *Name* and a *Description*. After creating a new Instance you can set it as current via the index page by clicking on the *Edit → Set current* button

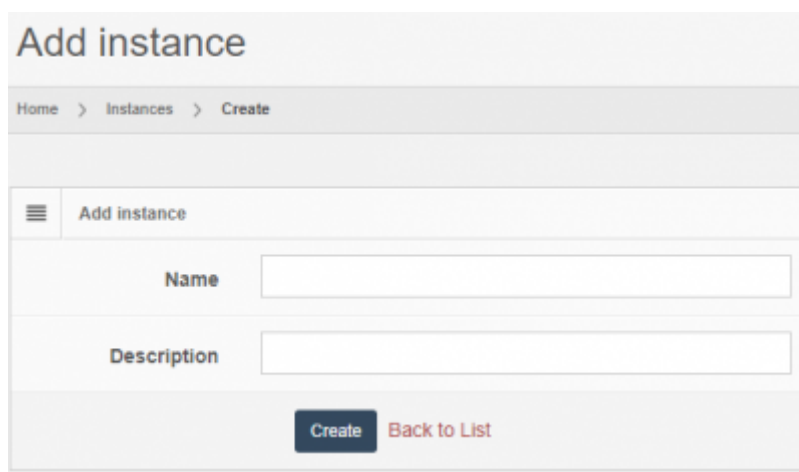

Add instance

Home > Instances > Create

Add instance

Name

Description

Create Back to List

From:

<https://docs.pathomation.com/pma.core/2.0.0/> - **PMA.core 2.x**

Permanent link:

https://docs.pathomation.com/pma.core/2.0.0/doku.php?id=settings_instances

Last update: **2022/04/01 18:03**

