

External script

External scripts is the functionality provided by PMA.core to configure and execute command line scripts via PMA.core itself.




Adding a new script

To configure a new script you need to provide the following information describing the script you want to execute and the parameters passed to it

Setting	Description
Name	A unique name for the script used to fetch it and distinguish from other
Command	The command line to execute, a bat file, cmd file or exe
Arguments	The arguments passed to the script
Parameters	A dynamic list of parameters passed to the script by PMA.core

The following types of parameters are supported

Parameter	Description
Number	A number parameter
Path	A path parameter i.e. a slide path
String	A string parameter

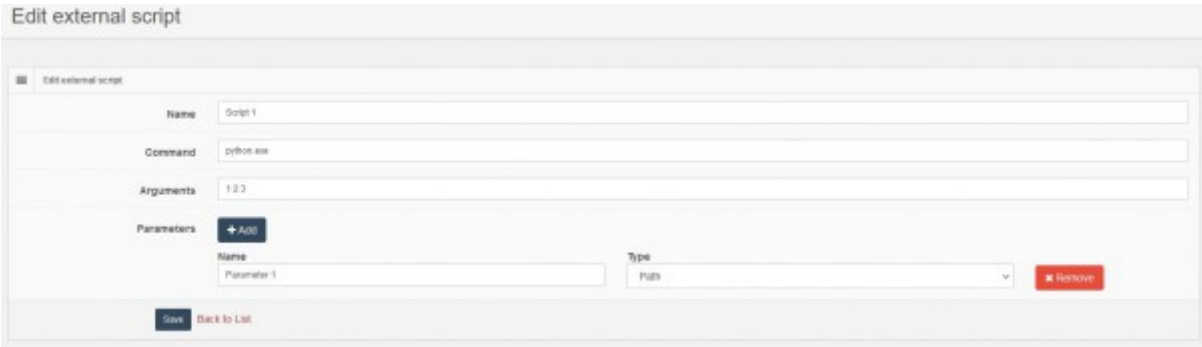

Executing an script

To execute an external script after adding and configuring it you have to utilize the [Scripts Run API](#)

Last update: 2022/04/01
18:30

settings_scripts https://docs.pathomation.com/pma.core/2.0.2/doku.php?id=settings_scripts&rev=1648827005

From:

<https://docs.pathomation.com/pma.core/2.0.2/> - **PMA.core 2.x**

Permanent link:

https://docs.pathomation.com/pma.core/2.0.2/doku.php?id=settings_scripts&rev=1648827005

Last update: **2022/04/01 18:30**

