

Simple shapes


Below are the set of simple shapes that can be added to a annotation.



- Point
- Circle
- Line
- Arrow
- Ellipse
- Rectangle
- Text
- Landmark 1
- Landmark 2
- Landmark 3


Point Shape

Lets start with drawing a point, once you select the shape and click on the slide; a point will be drawn and a popup dialog would appear that would allow you to add annotation title and classification. Once you press "Add Annotation" the annotation will be saved and will appear on the slide.

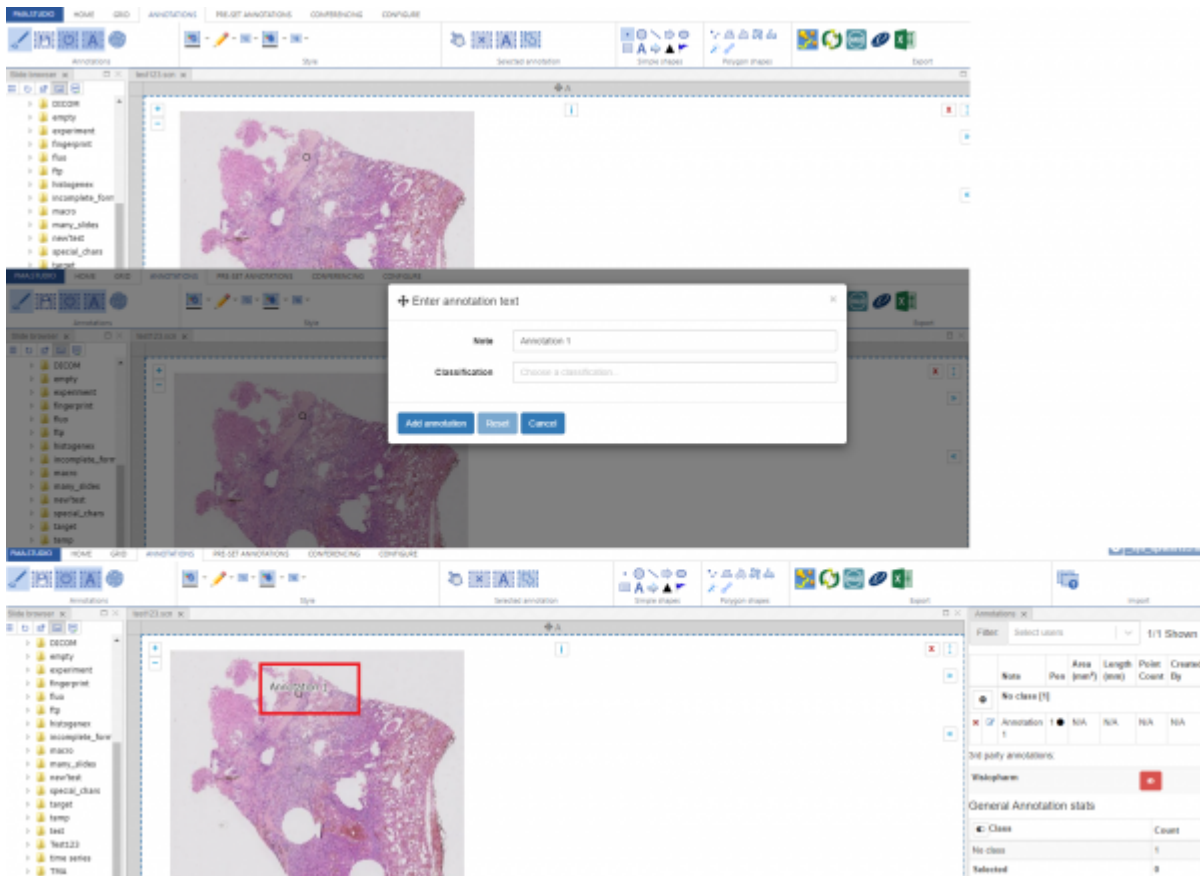


Circle Shape

Next is “Circle” tool that allows you to draw circle annotation on a slide.

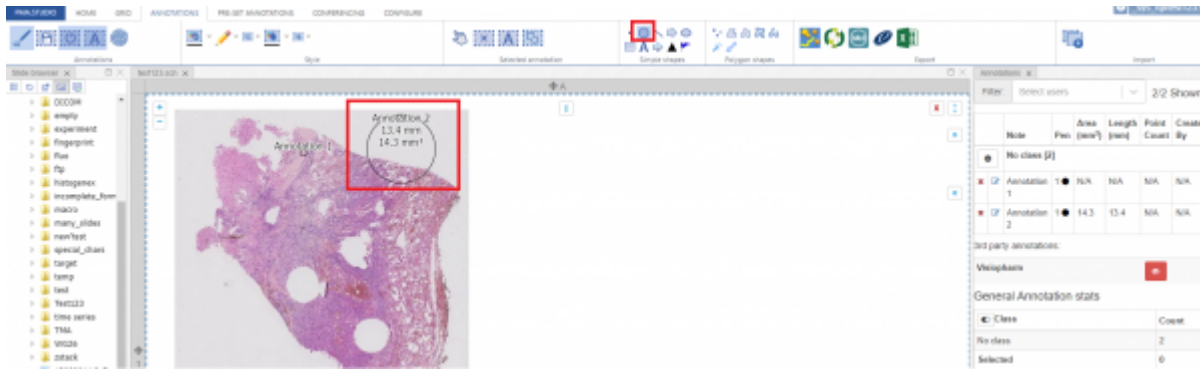


Line Tool

Next is “Line” tool that allows you to draw line annotation on a slide.

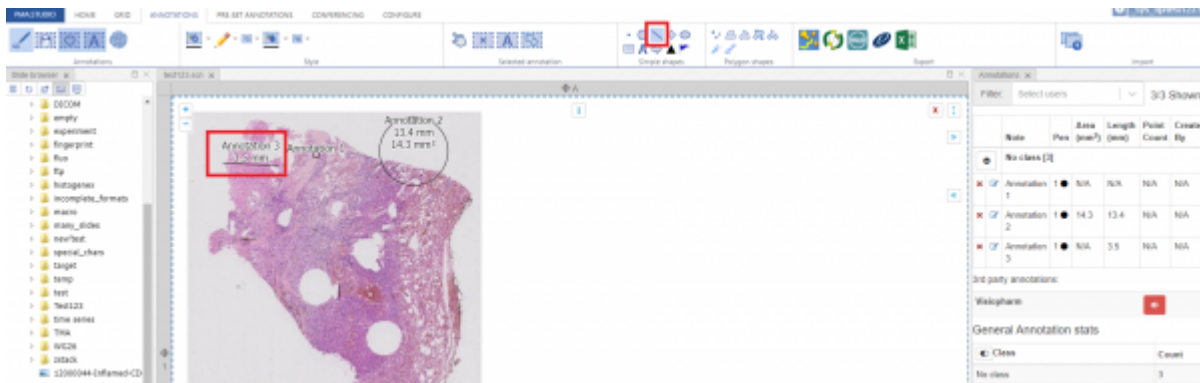


Arrow Tool

Next is “Arrow” tool that allows you to draw arrow annotation on a slide.

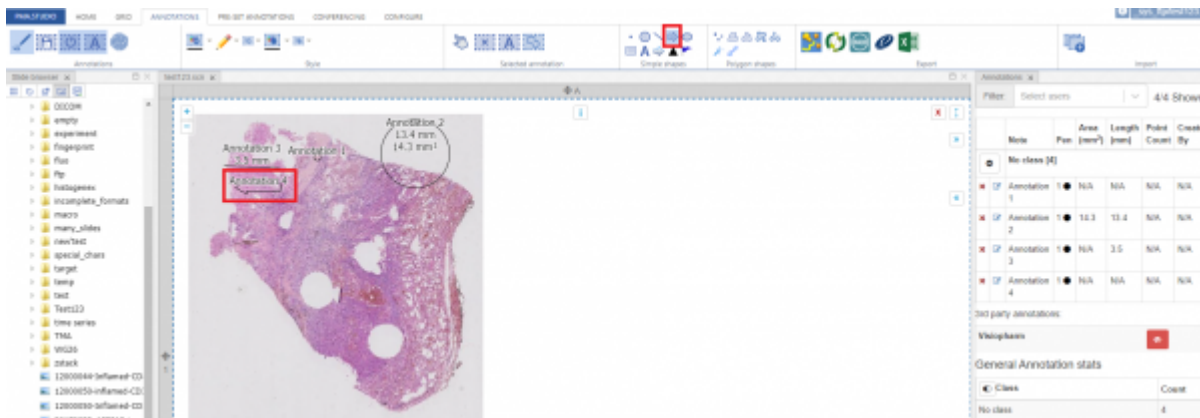

Ellipse Tool

Next is “Ellipse ” tool that allows you to draw ellipse annotation on a slide.

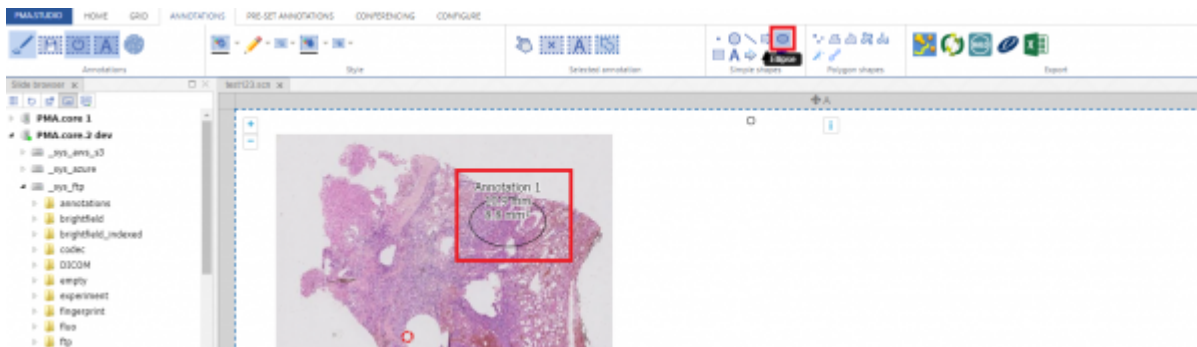


Rectangular Tool

Next is “Rectangular” tool that allows you to draw rectangular annotation on a slide.




Landmark 1 Tool

Next is “Landmark 1” tool that allows you to draw first of the 3 landmark annotation on a slide.

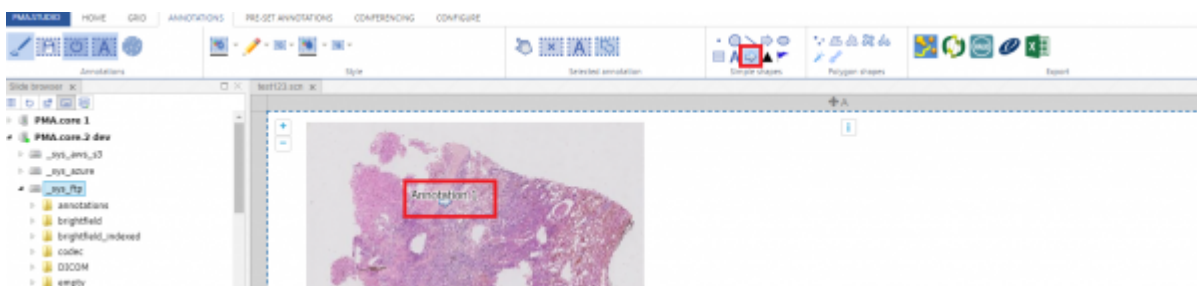


Landmark 2 Tool

Next is “Landmark 2” tool that allows you to draw 2nd of the 3 landmark annotation on a slide.

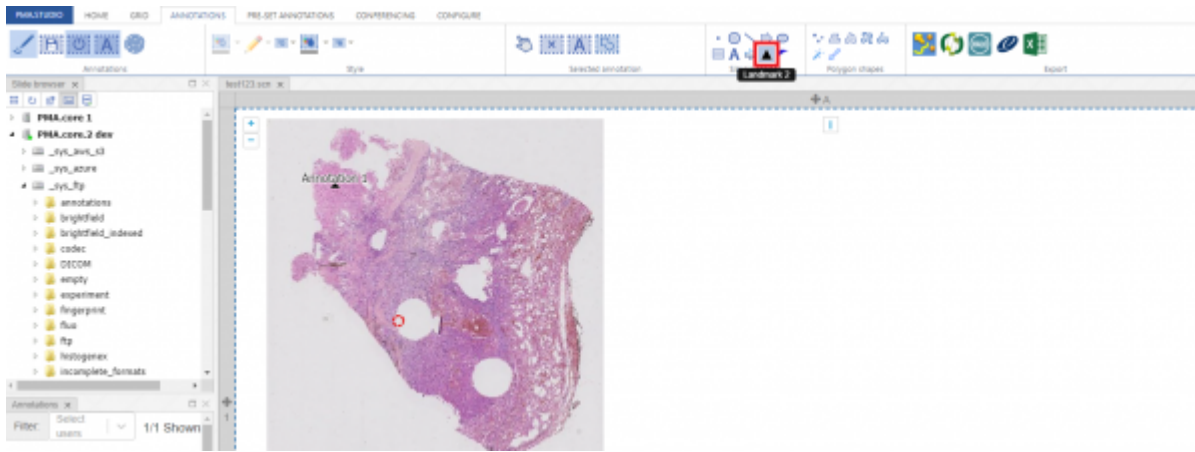

Landmark 3 Tool

Next is "Landmark 3" tool that allows you to draw last of the 3 landmark annotation on a slide.

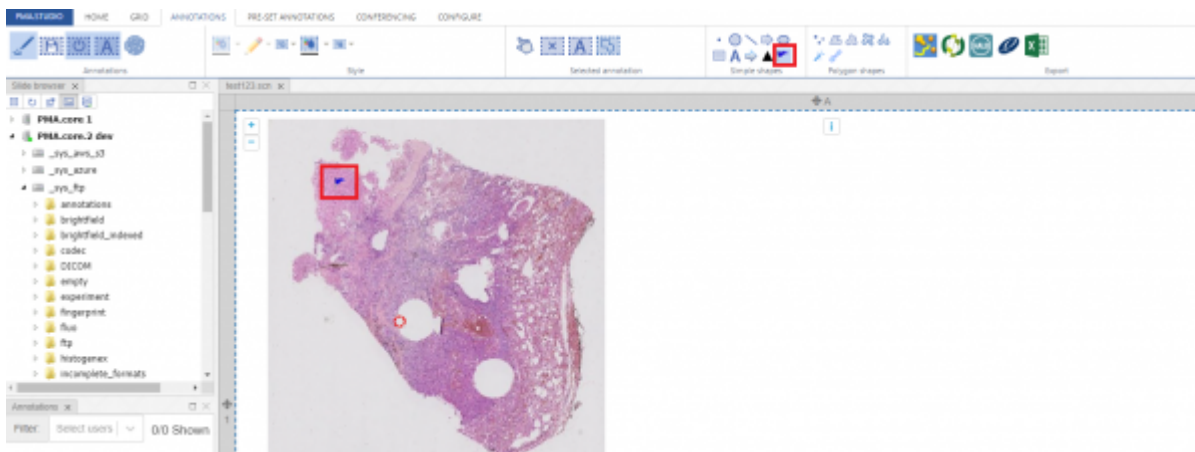

From: <https://docs.pathomation.com/pma.studio/2.1.0/> - **PMA.studio 2.1.0**

Permanent link: https://docs.pathomation.com/pma.studio/2.1.0/doku.php?id=annotations_simple&rev=1650555309

Last update: **2022/04/21 18:35**

