

## Forms panel

The Forms data panel purpose is to show filled forms for the slide and provide access to editing and filling new forms for currently selected slide. The panel consists of 2 main sections - Add button that will add a panel with specific form, and section with already filled form for the selected slide, user can collapse\expand each form with toggle control in top left corner and if user need to edit the form (s)he should click on 'Edit' button, it will open a new panel - specific form edit panel, where user can fill or edit form values and save them.

From:

<https://docs.pathomation.com/pma.studio/2.1.0/> - **PMA.studio 2.1.0**

Permanent link:

[https://docs.pathomation.com/pma.studio/2.1.0/doku.php?id=panel\\_forms&rev=1677772896](https://docs.pathomation.com/pma.studio/2.1.0/doku.php?id=panel_forms&rev=1677772896)

Last update: **2023/03/02 19:01**

